



# New Sterileware Sampling Tools


## Now Available!



Tools for Sampling from Bel-Art – SP Scienceware

### Innovative, Disposable Sterile Tools for Efficient and Reliable Sampling

- Reduce the risk of human error and sample contamination
- Made from FDA and EU conforming materials; safe for food, drug and cosmetic use
- Gamma irradiated sterile (SAL  $10^{-6}$ )
- Lot stamped for accurate tracing; shipped with a Certificate of Processing for Sterilization

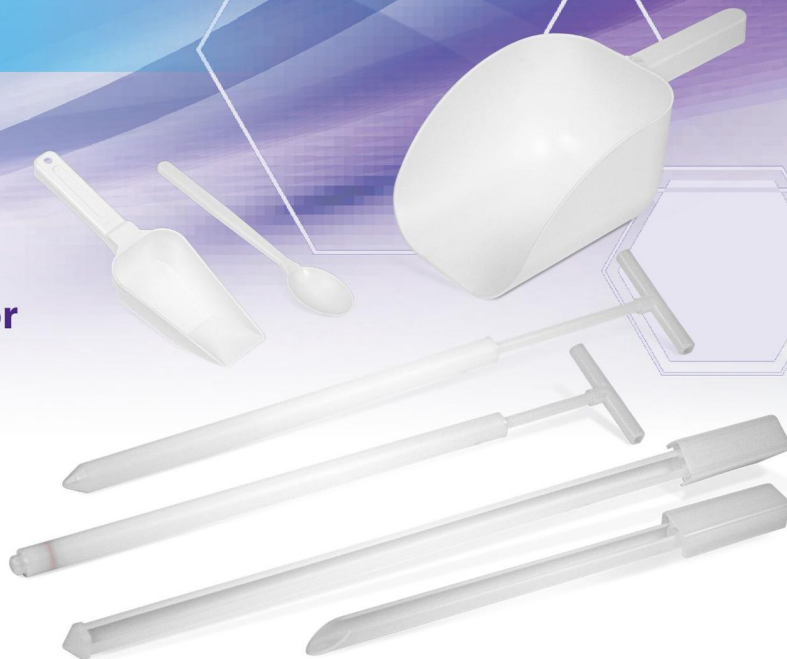


#### Sterileware Liquid Thief

##### Easily Sample Low Viscosity Liquids

CATALOG NO.	CAPACITY	QUANTITY/PACK
F37871-0500	100ml (3.4oz)	20
F37871-1000	190ml (6.4oz)	20

##### How it works:

1. Insert the sampler into the product.
2. At the required depth pull up the handle. The sample will be drawn into the sampler.
3. Withdraw sampler.
4. Push down handle to expel the sample.

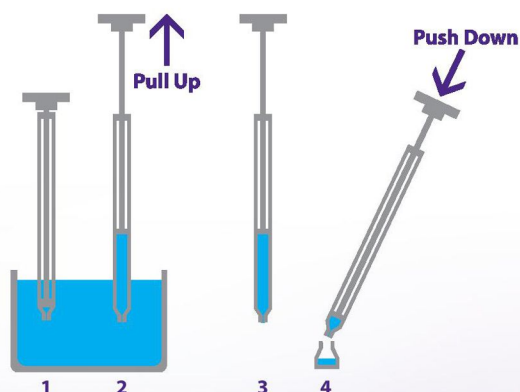


#### Sterileware Powder Thief

##### Sample from Within Bulk Powders

CATALOG NO.	CAPACITY	LENGTH	QUANTITY/PACK
F37874-0001	1ml (0.03oz)	50cm (19 <sup>5</sup> / <sub>8</sub> "	20
F37874-0005	5ml (0.17oz)	50cm (19 <sup>5</sup> / <sub>8</sub> "	20
F37874-0010	10ml (0.34oz)	50cm (19 <sup>5</sup> / <sub>8</sub> "	20
F37874-0020	20ml (0.67oz)	50cm (19 <sup>5</sup> / <sub>8</sub> "	20
F37874-1001	1ml (0.03oz)	100cm (39 <sup>3</sup> / <sub>8</sub> "	20
F37874-1005	5ml (0.17oz)	100cm (39 <sup>3</sup> / <sub>8</sub> "	20
F37874-1010	10ml (0.34oz)	100cm (39 <sup>3</sup> / <sub>8</sub> "	20
F37874-1020	20ml (0.67oz)	95.5cm (37 <sup>3</sup> / <sub>8</sub> "	20

##### How it works:

1. Insert the sampler into the product. Ensure that the tip is inside the sampler body.
2. At the required depth pull up the body to expose the tip. The sample will fall in around the tip.
3. Push down body of the sampler to trap the sample.
4. Withdraw sampler.
5. Pull up body to empty sample.

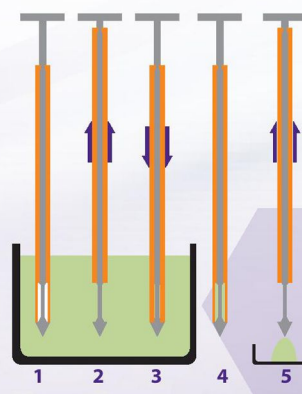

Available at



**DAWN  
SCIENTIFIC INC**  
LABORATORY SUPPLIES & EQUIPMENT



Tools for Sampling from Bel-Art – SP Scienceware



### Sterileware Pharma Scoops - White High Quality, Polystyrene Scoops



CATALOG NO.	NOMINAL CAPACITY	QUANTITY/PACK
F36909-0030	30ml (1oz)	100
F36909-0060	60ml (2oz)	100
F36909-0125	125ml (4oz)	100
F36909-0250	250ml (8oz)	85
F36909-0500	500ml (17oz)	50
F36909-1000	1000ml (34oz)	25
F36909-2500	2500ml (85oz)	15

### Sterileware Flat Bottom Sampling Scoops – White Useful as a Weigh Boat or for Volume Measurements



CATALOG NO.	NOMINAL CAPACITY	QUANTITY/PACK
F36911-0025	25ml (0.8oz)	100
F36911-0050	50ml (1.7oz)	100
F36911-0100	100ml (3.4oz)	100
F36911-0150	150ml (5.1oz)	100
F36911-0250	250ml (8oz)	100
F36911-0500	500ml (17oz)	40
F36911-1000	1000ml (34oz)	30

### Sterileware Oval Sampling Spoons – White Small Sample Collection; Long Handle Reach



CATALOG NO.	NOMINAL CAPACITY	QUANTITY/PACK
F36941-0003	3ml (0.1oz)	100
F36941-0010	10ml (0.3oz)	100

### Sterileware Powder Lance Samplers Cohesive Powder Sampling Solution



CATALOG NO.	CAPACITY	QUANTITY/PACK
F36974-0420	75ml (2.5oz)	20
F36974-0600	125ml (4oz)	20
F36974-0950	200ml (6.8oz)	20

### Sterileware Powder Probe Samplers Free-Flowing Powder Sampler



CATALOG NO.	CAPACITY	QUANTITY/PACK
F36975-0600	125ml (4oz)	20
F36975-1000	200ml (6.8oz)	20



WARNING – Important information for California residents regarding Prop 65. Visit [www.P65Warnings.ca.gov](http://www.P65Warnings.ca.gov) for more information.

Available at :



121 LIBERTY STREET, METUCHEN, NJ, 08840

TOLL-FREE NUMBER  
1-800-DAWN-SCI

FOR INQUIRY  
+1 732-902-6300  
SALES@DAWNSCIENTIFIC.COM  
WWW.DAWNSCIENTIFIC.COM

